

SLIT DISK REPLACEMENT

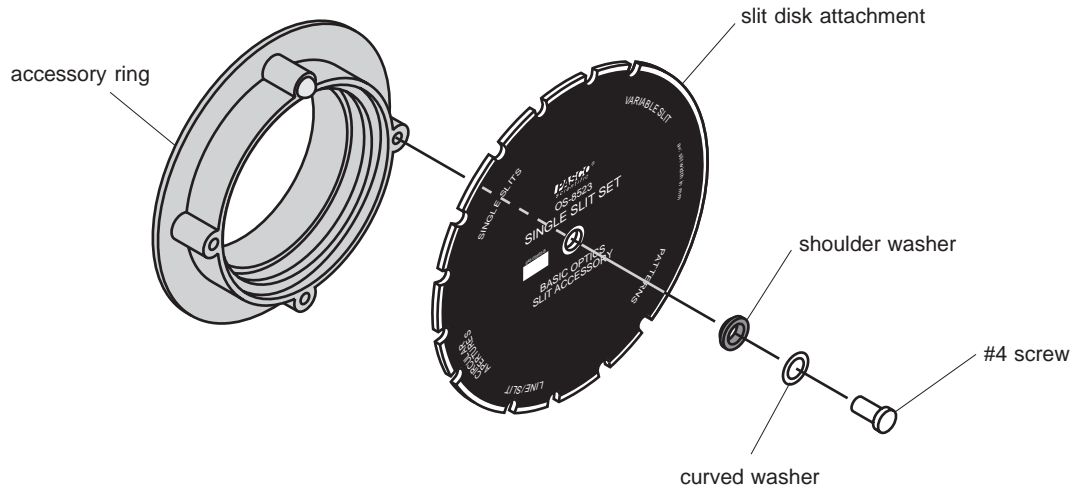


Figure 1: Slit Disk Attachment Replacement

Introduction

These instructions allow the user to replace Single and Multiple Slit disk attachment plates for the PASCO OS-8523 Slit Accessories.

Equipment

This kit includes the following parts:

- Single Slit disk
- Multiple Slit disk
- #4 X 3/8" self-tapping screws (2)
- 3.5mm (inside diameter) curved spring washers (2)
- 3mm (inside diameter) shoulder washers (2)

You will need the following tools:

- #1 Phillips-drive screwdriver

Assembly

- ① Remove old or damaged slit disk attachments. Please discard attachments.
- ② Assemble supplied replacement slit disk attachments as shown in Figure 1. Use supplied hardware if necessary. Slit disks must rotate yet still "lock" into position. Note orientation of accessory rings. **DO NOT OVERTIGHTEN SCREWS.**
- ③ Please refer to the instruction manual shipped with your OS-8523 Slit Accessories.

© 1996 PASCO scientific

This instruction sheet written/edited by:

Mary Ellen Niedzielski

Notes: